

Hygiene & Environmental Testing Pro-media **SWAB TEST**

(Patented)

Diluent

PBS

Phosphate Buffered Saline

BPW

Buffered Peptone Water

MRD

Maximum Recovery Diluent



SWAB TEST kit has been certified
by HACCP International.
(Asia, Oceania)



All in One

Ready to use self
contained swab and
diluent system.

No pipetting require

Flip the lid and inoculate media
using the measuring
gauge on the side.

Extension Tool

Effective to sampling
for hard to reach
places.

No test tube require

Incubate the entire
container after
added a media
ST-SSP for coliform
group test.

Compact incubator allows swabs to be readily processed on-site.

Manufacturer **ELMEX Limited**

2-2 Ichigayasadohara-cho, Shinjuku-ku, Tokyo 162-0842 JAPAN
TEL : 081-3-5229-7911 FAX : 081-3-5261-4130
URL : <http://www.elmex.co.jp> e-mail address : sales@elmex.co.jp

Basic Operation



Squeeze swabbud to remove excess diluent.



Swab.



Mix to suspend.



Squeeze to dispense.

Suggested Use

<Ex.1> Quantitative test

Flip open the lid, invert and inoculate by squeezing onto your media.

Each Swab test contains 10mL of diluent and the dispensed volume can be measured using the gauge on the side of the bottle.

Alternatively, one drop is 50µL or 20 drops per mL. The screw cap can be removed to allow the pipetting directly from the tube.




<Ex.2> Qualitative test (Presence or Absence)

For Coliform group test, simply add the ST-SSP reagent directly into Swab Test tube via the dispensing nozzle. Incubate for 24 hours at 36°C.



<Result>

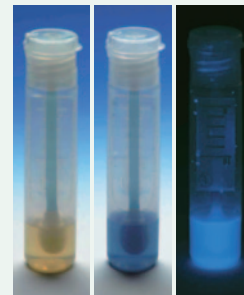

Coliform G.	-	+	+
<i>E.coli</i>	-	-	+

※ Refer to its instruction about culture media

Extension Tool

(optional)


Use an extension tool to swab hard-to-reach areas.

Insert the stick into the capper and attach it to the cap of the SWAB TEST kit before use. Stick is able to concatenate 1, 2 or 3 pieces.



capper

stick

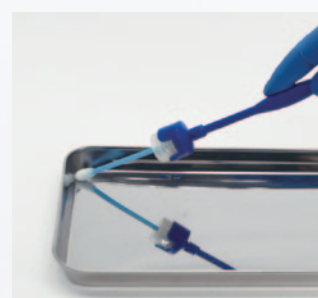

Code No.	Quantity/box	Length (mm)	Material
SE-4	4 pairs (capper × 4pcs) stick × 4pcs)	Capper + Stick × 1pcs = 155 Capper + Stick × 2pcs = 280 Capper + Stick × 3pcs = 405	ABS

Code No.	ST-25PBS	ST-25BPW	ST-25MRD
Diluent	Phosphate Buffered Saline (0.85% NaCl)	Buffered Peptone Water (1.0% Peptone)	Maximum Recovery Diluent (0.1% Peptone)
Quantity/box	400pcs (10pcs×40)		
Storage	Room temperature (2~25°C) , protected from direct light		
Shelf life (Unpacked state)	Two (2) years from production date.		
Sterilization	Gamma irradiation		

◆ Specifications are subject to be changed for enhancement, with out notice.

Manufacturer **ELMEX Limited**

2-2 Ichigayasadohara-cho, Shinjuku-ku, Tokyo 162-0842 JAPAN
TEL : 081-3-5229-7911 FAX : 081-3-5261-4130
URL : <http://www.elmex.co.jp> e-mail address : sales@elmex.co.jp